

New and Unique Techniques in Life Science using Microplate Instrumentation from BioTek Instruments

Different Cell-Based Assays using the Patented Hybrid Technology™ and Cell Imaging

Dr. Andreas Rieger¹,

¹BioTek Instruments, Inc., Winooski, VT, USA

BioTek Instruments has been awarded US patent 8,218,141 covering the combination of two alternate paths for light within the same analytical instrument. Known as Hybrid Technology™, the patented functionality is available in BioTek's Synergy™ NEO HTS Multi-Mode Microplate Reader and in the Synergy H4 and H1 Hybrid Multi-Mode Microplate Readers.

Using Hybrid Technology™, light can be directed to and from the microplate samples through a filter-based optical system or through a monochromator-based optical system for wavelength selection. This two-optical-path system provides optimal performance and flexibility for a wide range of applications in fluorescence microplate assays. BioTek's Synergy readers are modular platforms that can include Fluorescence Intensity, UV-Vis Absorbance, Luminescence, TRF, TR-FRET and FP detection modes.

This presentation we will demonstrate the power of BioTek instrumentation which can perform a wide range of diverse applications in the biochemistry laboratory. These applications cover micro-volume nucleic acid quantification from 2 µL samples using the monochromators in absorbance mode; high sensitivity cellular signal transduction assays (AlphaScreen™) using spectral filters; live cell fluorescence assays using the monochromators in fluorescence mode and finally using the GloSensor® (Promega) for monitoring cAMP levels of GPCRs using a bioluminescent readouts or powerful Imaging to be used for phenotypic analysis or cell counting.

Hybrid Technology™

Combined systems



▪ **Combined benefits of filter-based and monochromator-based systems:**

- Flexibility and convenience of monochromator-based systems
- Sensitivity and speed of filter-based systems

